**Xavier Garcia Elena**

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Balmes, 195, 7º 2ª - 08006 Barcelona  
Tel. 93 415 99 90 - info@ampersand.net

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**Certification**

Mr. Xavier Garcia Elena, Official Translator of the English, French and Catalan languages, certifies that the translation contained in the previous pages is a true and complete rendering into Spanish of a document written in English.

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Cells4Life Group LLP – Form 100 - SOP 100 Template v6			
SOP Name	Receipt of Samples		
SOP Number	112		
SOP Version No.	v. 5	Supersedes Version No.	v. 3
SOP Issue Date	August 2012	SOP Revision Date	August 2013
SOP Written by (initials)	RED	SOP Authorised by (initials)	RR
<b>DO NOT copy, amend or distribute without prior written authorisation from Cells4Life</b>			

Scope of SOP	The aim of this SOP is to describe the correct action on receipt of UCB and UCT samples at the Cells4Life laboratory
Other Documents Referred to	Sugar Operating Manual
Risk Assessments Referred to	

1. Samples will be delivered to the Cells4Life laboratory by both local and international courier under appropriate agreement with Cells4Life.
2. When a sample is delivered proof of delivery must accompany the sample.
3. When a kit is delivered to the Cells4life the kit is to be taken directly to laboratory reception.
4. Inspect sample boxes/packages for obvious signs of damage and/or hazards before handling..
5. If any such hazards are observed immediately notify a member of lab staff and stop other people from handling the sample.
6. Persons signing for the sample MUST enter the kit number and the time delivered to laboratory is to be noted in the kit receipt diary under the date of reception, (found on reception) and enter the following:
  - i. Sample/Kit number – the unique sample identifier
  - ii. TTL (time to lab) – the time the sample was signed for
  - iii. TP (time processed) – for lab staff to complete
7. All UK-based samples must be placed in the fridge in lab reception
8. All non-UK samples i.e. ‘NanoCool’ boxes can be left on the bench in reception
9. If laboratory staff is present, notify the member of laboratory staff that the sample has arrived.
10. If lab staff are not present email ‘courier@cells4life.co.uk’ or the individual lab person on-call to notify them of the sample arrival.
11. All associated paperwork must be left in the appropriate tray in lab reception.

Note:- The following steps are to be carried out only by trained laboratory staff.

12. Ensure you are wearing a laboratory coat, goggles and gloves.
13. Always handle all sample as if contaminated, even if its status is known.
14. If the kit has been transported in an outer transportation box, remove it from the transportation box.
15. Switch the webcam to record.
16. Record the specimen box, noting if the box is sealed or not, ensure that the kit number is in focus.
17. Open the box.
18. Ensuring that there are no sharps in the box, (i.e. needles), carefully remove the contents of the box individually showing each item to the webcam. Ensure that the kit number is in focus. If any neeles are present remove and place in sharps container.
19. Confirm that the box is empty by showing the inside to the webcam.
20. Stop recording.

Cells4Life Group LLP – Form 100 - SOP 100 Template v6			
SOP Name	Receipt of Samples		
SOP Number	112		
SOP Version No.	v. 5	Supersedes Version No.	v. 3
SOP Issue Date	August 2012	SOP Revision Date	August 2013
SOP Written by (initials)	RED	SOP Authorised by (initials)	RR
<b>DO NOT copy, amend or distribute without prior written authorisation from Cells4Life</b>			

21. Rename the video file to the corresponding kit number.
22. Place the specimen(s) into a specimen tray, (one kit number per tray), and place into the specimen reception fridge, until you are ready to process the sample(s).
23. Log the sample onto Sugar – refer to Sugar Operating manual.
24. Convert the video file from an .msv to an .avi file.
25. Upload the unpacking video to Sugar, attached to the birth record. Refer to the Sugar Operating Manual.
26. Any issue associated with the kit including the manner in which it was packed is to be noted in Sugar under kit birth record.